

USER INSTRUCTIONS

DRYING CABINETS



THIS APPARATUS MUST BE EARTHED.

For a complete electrical disconnection pull out the mains plug.

On arrival unpack carefully.

Place cabinet on an even, level surface ensuring that the door(s) are level.

Fit the shelves at the required positions.

With sliding door models put glass into top runner and then into lower runner.

Plug the cabinet into a suitable power outlet.

Turn the control dial to the appropriate heating power setting.

The green power lamp and amber heat lamp will illuminate.

The amber lamp will switch on and off as the elements are controlled.

Ensure that the shelves are not blocked with trays, paper materials, etc.

THIS CABINET IS UNSUITABLE FOR DRYING ITEMS THAT HAVE BEEN RINSED WITH ANY TYPE OF ALCOHOL OR SOLVENT.

Cleaning.

This is an electrical appliance so the electrical supply must be disconnected first.

The chamber should only be cleaned with a damp cloth.

If you have any concerns or want additional advice please contact your supplier: